


INSTRUCTIONS FOR ASSEMBLY

MODEL C-10 SERIES CHROMATO-VUE® CABINET

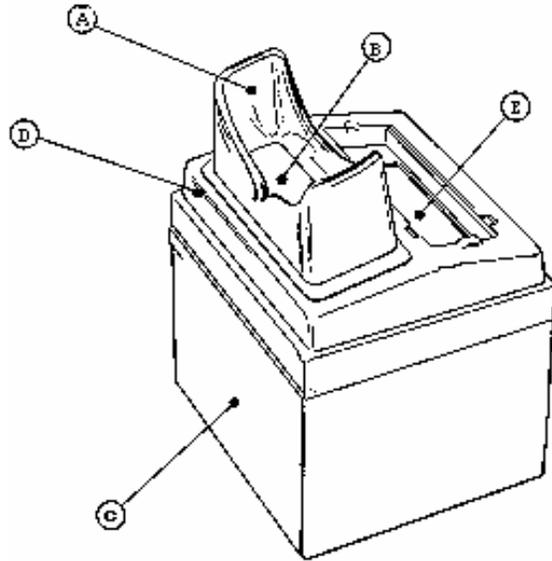

1. Lift flexible eyepiece (A) and cover assembly (B) out through top of cabinet.
2. Slip lamp well (E) end of cover assembly over rim at back of cabinet until flush with top. Apply enough pressure towards front portion to snap over rim. Check to see that cover assembly is flush with top on all sides.
3. Slip flexible eyepiece over rim above contrast control filter (D) with notched portion facing front of cabinet.
4. Place your Mineralight lamp in molded lamp well with line cord at left side. Your Chromato-Vue cabinet is now ready to use. Lift light curtain (C) when placing or positioning materials; leave in closed position when in use to block out any external light.

 UVP

C-10P

CHROMATO-VUE CABINET

The difference between the C-10, C-10E and CC-10P Cabinets is the addition of a lamp support bracket on the C-10P. The purpose of the bracket is to hold the portable lamp securely in place.

To assemble the C-10P, follow the C-10 Cabinet Series instructions numbers 1 through 3. Instruction No. 4 for the C-10P follows:

4. Place your portable lamp in the lamp support bracket. This bracket firmly holds the lamp in place. Your Chromato-Vue cabinet is now ready to use. Lift light curtain (C) when placing or positioning materials; leave curtain in closed position when in use to block out any outside light.

Replacement parts Cabinets:

Eyepiece: 23-0036-01
Contrast Control Filter: 38-0013-01



Corporate Headquarters: UVP, LLC

2066 W. 11th Street Upland, CA 91786 / Tel: (800) 452-6788
(909) 946-3197 / Fax: (909) 946-3597

European Operations: Ultra-Violet Products Ltd

Unit 1, Trinity Hall Farm Estate, Nuffield Road,
Cambridge CB4 1TG UK / Tel: +44(0)1223 420022
Fax: +44(0)1223 420156

For additional information on UVP's products, go to www.UVP.com

81-0008-01 Rev F